

## Quaternary amines analysis by LCM and finding the right ion -How To

When analyzing quaternary amines in LC-MS, you should remember that they are permanently charged species.

A tertiary amine for example, may be observed in the positive ion mode as the [M + H]<sup>+</sup> ion, but because quaternary amines are already ionized, you should look for the [M]<sup>+</sup> ion in the extracted ion chromatograms (EICs) instead.

If you look for the m/z of the  $[M + H]^+$  ion, you may see no peak or a peak corresponding to some other compound in the sample because quaternary amines do not form these ions.

These kinds of suggestions can help you identify the proper peak for your <u>analyte</u> and avoid peak assignment errors.



Printed from the Chrom Resource Center Copyright 2025, All Rights Apply **MicroSolv Technology Corporation** 9158 Industrial Blvd. NE, Leland, NC 28451 Tel: (732) 380-8900 Fax: (910) 769-9435 Email: customers@mtc-usa.com

Website: www.mtc-usa.com