



Since Cogent TYPE-C columns can be used in Reversed Phase or Normal Phase what do you suggest to switch these solvent systems which are not miscible - FAQ

Suggested Procedure:

A – Changing from Reversed Phase to Normal Phase HPLC; pump 100% methanol for 15 minutes at 1 mL / minute flow rate, followed by 15 minutes 100% methylene chloride. The column is ready to be equilibrated with mobile phase for NP HPLC.

B – Changing from Normal Phase to Reverses Phase HPLC; pump 100% methylene chloride for 15 minutes at 1 mL / minute flow rate, followed by 15 minutes 100% methanol. The column is ready to be equilibrated with mobile phase for RP-HPLC.

A simple 30 minute procedure allows switching from one mode to another.



Printed from the Chrom Resource Center
Copyright 2025, All Rights Apply
MicroSolv Technology Corporation
9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900
Fax: (910) 769-9435
Email: customers@mtc-usa.com
Website: www.mtc-usa.com