

Can autosampler vials be used in a centrifuge and what is the maximum centrifugal force that can be used - FAQ

Autosampler vials are not designed to be centrifuged so we cannot support it with the exception of the MRQ™ & Max Recovery™ vials.

Due to their thicker walls and "single piece" construction (AQ brand, RSA or RSA-Pro) they can withstand g forces up to 6,000 x g when used in a compatible centrifuge and adapters.

Click [HERE](#) for ordering information and pictures of the MRQ™ center draining vials.

Click [HERE](#) for ordering information and pictures of the Max Recovery™ vials.

Click [HERE](#) for ordering information and pictures of The Vial Centrifuge product page

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com