


## Hydrophobic and Hydrophilic Peptides Analyzed with HPLC - AppNote

### Simultaneous Separation using a Capillary Column

The studied analytes comprised some very Hydrophobic and Hydrophilic Peptides with diverse isoelectric points (pI) and thus served as useful probes to explore the potential of this App Note in this field of Bioanalysis.

The figure shows how these Peptides of widely disparate physicochemical properties could be separated in a single run. With these capabilities, laboratories can be more efficient and increase throughput by not having to perform separate analyses for both Hydrophobic and Hydrophilic Peptides using different Columns.

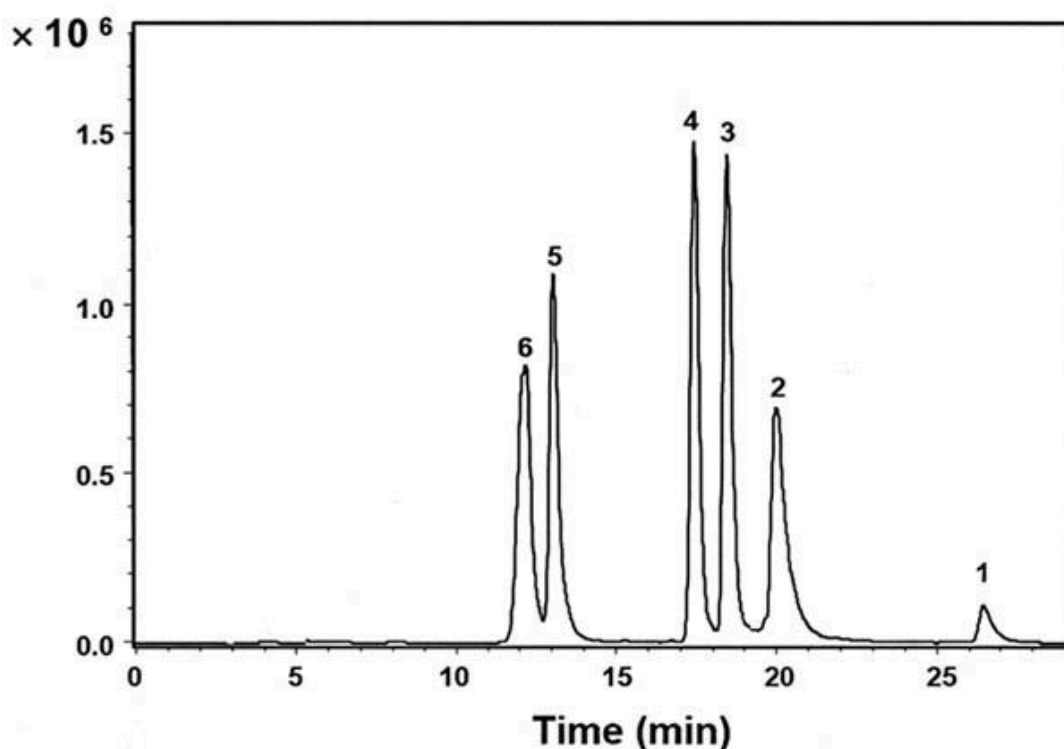

Peptide code	Sequence	Calculated hydrophobicity	pI
1	H-Gly-Arg-Ala-Asp-Ser-Pro-Lys-OH	-2.31	9.4
2	H-Lys-Gln-Ala-Gly-Asp-Val-OH	0.35	5.8
3	H-Tyr-Ile-Gly-Ser-Arg-OH	4.11	9.1
4	H-Trp-His-Trp-Leu-Gln-Leu-OH	9.65	7.1
5	H-Tyr-Gly-Gly-Phe-Leu-OH	10.61	5.5
6	H-Tyr-Tyr-Tyr-Tyr-Tyr-Tyr-OH	11.34	5.5

### Peaks:

1. H-Gly-Arg-Ala-Asp-Ser-Pro-Lys-OH
2. H-Lys-Gln-Ala-Gly-Asp-Val-OH
3. H-Tyr-Ile-Gly-Ser-Arg-OH
4. H-Trp-His-Trp-Leu-Gln-Leu-OH
5. H-Tyr-Gly-Gly-Phe-Leu-OH
6. H-Tyr-Tyr-Tyr-Tyr-Tyr-Tyr-OH

### Method Conditions:

**Column:** Cogent Diamond Hydride™, 4um, 100A R&D Capillary Column

**Dimensions:** 0.3mm x 150mm

#### Solvents:

A: DI Water / 0.5% Formic Acid (v/v)

B: Acetonitrile / 0.5% Formic Acid (v/v)

#### Gradient:

Time (minutes)	%B
0	90
5	90
10	70
20	60
20.1	30
30	30

**Injection Volume:** 0.1µL

**Flow Rate:** 4µL / minute

**Detection:** ESI - POS - Agilent 6210 MSD TOF Mass Spectrometer

**Sample Preparation:** 10µg / mL of each Peptide in 50% Solvent A / 50% Solvent B.

**Notes:** The data and discussion presented here are delineated in greater detail in a full research article. Reference: R.I. Boysen, Y. Yang, J. Chowdhury, M.T. Matyska, J.J. Pesek, M.T.W. Hearn, "Simultaneous separation of hydrophobic and hydrophilic peptides with a silica hydride stationary phase using aqueous normal phase conditions." *J. Chromatogr. A*, 2011, 1218, 8021–8026.



### Attachment

No 371 Hydrophobic and Hydrophilic Peptides Analyzed with HPLC pdf 0.3 Mb [Download File](#)

Printed from the Chrom Resource Center  
Copyright 2025, All Rights Apply  
**MicroSolv Technology Corporation**  
9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900  
Fax: (910) 769-9435  
Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)  
Website: [www.mtc-usa.com](http://www.mtc-usa.com)