


## Glutathione (GSH) Analyzed by LCMS - AppNote

## **LCMS Method without Derivatization**

The major benefit of this analysis method for Glutathione which uses the Cogent Diamond Hydride Column and MS detection is that it is performed without derivatization, unlike other LCMS approaches. A GSH signal of MS is stable after repeated analyses. The unique design of the ESI offers efficient supersoft ionization which helps to preserve the structure of the analyte.

This Method ensures low variability of the results. After validation, it could be applied to routine analysis of oxidative stress in clinical samples (blood or plasma).

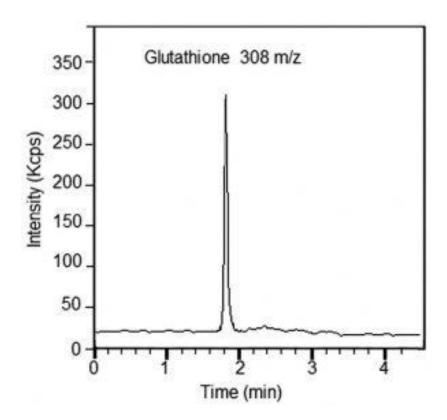

Glutathione

Peak:

Glutathione 308 m/z

## **Method Conditions**

Column: Cogent Diamond Hydride 2.o™, 2.2µm, 120Å

Catalog No.: 70200-05P-2

Dimensions: 2.1 x 50mm

**Mobile Phase:** 

A: DI Water / 0.1% Formic Acid (v/v)
B: Acetonitrile / 0.1% Formic Acid (v/v)

**Gradient:** 

Time (minutes)	%B
0	85
2	30
4	30
5	85

**Post Time:** 3 minutes **Flow rate:** 0.3 mL/minute

**Detection:** ESI – POS - PerkinElmer Flexar SQ 300 Mass Spectrometer

Injection vol.: 1µL

Sample Preparation: 1 ppm of Glutathione in 50% Acetonitrile / 50% DI Water / 0.1% Formic Acid

**t**<sub>0</sub>: 0.7 minutes

Note: Glutathione (GSH, Gamma-L-Glutamyl- L-Cysteinyl-Glycine) is a non-protein Thiol compound which plays an important role in protecting calls against oxidative stress. Lower levels of Glutathione were found in different diseases, like cardiovascular disease, cancers, and aging [3,3]. Under oxidative stress, GSH is converted into its oxidized form Glutathione Disulfide (GSSG). A lower level of GSH is an early indicator of disease risk.

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com



## **Attachment**

No 279 Glutathione pdf 0.2 Mb Download File

Printed from the Chrom Resource Center
Copyright 2025, All Rights Apply
MicroSolv Technology Corporation
9158 Industrial Blvd. NE, Leland, NC 28451