

Is the Cogent UDC Cholesterol HPLC column stable as a stationary phase - FAQ

While cholesterol is a large molecule, the UDC-Cholesterol phase has proven to be stable in 0.1% v/v TFA and 0.1% v/v H₃PO₄ at acid pH's and up through pH 8.00 with 20mM NH₄COOH buffer with NH₄OH.

A pH in excess of pH 8.00 should be avoided as this phase has a mono dentate (*single point*) attachment to the silica matrix and an intermediate ester link in its structure.

For higher pH's the Cogent Bidentate C18™ is recommended.



Printed from the Chrom Resource Center
Copyright 2025, All Rights Apply
MicroSolv Technology Corporation
9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900
Fax: (910) 769-9435
Email: customers@mtc-usa.com
Website: www.mtc-usa.com